



MP Biomedicals, LLC

29525 Fountain Parkway
Solon, Ohio 44139

Telephone: 440/337-1200
Toll Free: 800/854-0530
Fax: 440/337-1180
mailto: biotech@mpbio.com
web: <http://www.mpbio.com>

TECHNICAL INFORMATION

Catalog Number: 821682, 821683, 821684, 821890
RNase Erase

Instructions For Use

RNase Erase™ was formulated using three ingredients known to be active against RNase. It effectively removes high levels of RNase contamination that similar products cannot. Additionally, RNase Erase™ has been formulated so that it can be used to remove RNase contamination from reaction vessels. With thorough rinsing, RNase Erase™ leaves no residues that are inhibitory to enzymatic reactions. The solution is ready to use. If there is a precipitate (as may happen at low temperatures), shake and/or heat at 37°C to bring the precipitate back into solution. The product should be stored at room temperature. Always wear gloves while using RNase Erase™ as prolonged contact with skin may cause irritation. RNase Erase™ is not compatible with corrodible metal surfaces.

Instructions

Cleaning work surface: Apply directly to surface to be cleaned, wipe thoroughly with paper towel, rinse with water and then dry with a clean paper towel.

Cleaning lab apparatus: Apply RNase Erase™ liberally to a paper towel and wipe all exposed surfaces of the apparatus thoroughly. Rinse with water and then wipe dry. Some small parts may be cleaned by briefly soaking them in RNase Erase™, rinsing them in water and then drying. Please note that RNase Erase™ should not be used on corrodible metal surfaces.

Cleaning plastic and glass vessels: Add enough RNase Erase™, so that the entire surface of the vessel can be coated with the solution upon swirling, or vortexing in the case of centrifuge and microfuge tubes. After discarding the solution, rinse vessels thoroughly two times with distilled water.

Cleaning pipetters: Remove shaft from pipetter according to the manufacturers instructions. Remove seals and gaskets from the shaft and then soak shaft for one minute in RNase Erase Rinse the shaft thoroughly with water and then reassemble pipetter

Warning: If contact with eyes occurs, immediately flush with water and call a physician. If ingested, do not induce vomiting, give plenty of water and contact a physician.

ORDERING INFORMATION:

Catalog Number	Description	Size
821682	Spray Bottle	250 ml
821683	Dropper Bottle	2 x 125 ml
821684	Refill Bottle	250 ml
821890	Spray Bottle/Refill	1 liter